Product Datasheet: Precision Series Piston Pumps

The Precision series piston pumps were developed specifically for applications requiring a very exact flow at a fast dispense rate. The piston geometry isolates the outlet and inlets without the need for internal check valves, simplifying the construction and reducing flow restriction. The pumps are typically provided in a "fixed" position to deliver a set flow-per-stroke. An adjustable mounting frame is optional, allowing the pump displacement to be altered in the field and then locked in position. Speed sensing options include encoders, hall sensors, and/or photo detectors. The displacement algorithm is well-defined, allowing the end user to correlate motor position with the amount of fluid dispensed.

All pumps are customized to suit your requirements: the information below represents one of the possibilities...



9mm Piston Pump with Adjustable Displacement

19mm Piston Pump with Adjustable Displacement

- Simple construction reduces cost
- Positive Displacement: accurate and repeatable flow rates
- Mounting Versatility: Mounting brackets are customized for each application
- Chemically inert wetted materials
- Select from five piston diameters
- Available with stepper, brush-less DC, and brush-type DC motors
- 100% tested to customer requirements






6.34mm Piston Pump with Electronic Dosing Control

QF3222i Product Datasheet



Piston pumps undergoing long-term life testing

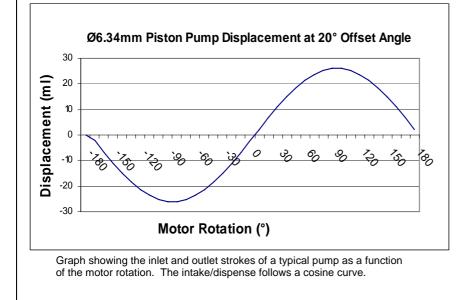

Specifications

Performance:

Max. continuous pressure: 4 bar Max. intermittent pressure: 6 bar Max. static case pressure: 6 bar Inlet: Self-priming Accuracy: within 1% of full scale Precision: Typically < 0.5% Speed range: 0-1500 rpm Fluid viscosity range: 0.3 to 1000 cps

Electrical:

Motor type: typically stepper motor, but also available in brush and brush-less DC. Rotation: Reversible Max. current: Dependent on motor type Max. output power: Dependent on motor type Connection: Flying wires, cable, or tabs Motor rating: Motor specific Speed sensing: Optional

Temperature: Fluid temp. range: -20 to +100°C Ambient air temp. range: 0-60°C Relative humidity range: 0-95% noncondensing

Construction: Standard: PVDF, Ceramic (alumina), PTFE. Optional: PEEK[®], Polysulfone, Polypropylene Rinse ports: Optional Inlet/Outlet: 3mm barbed tube (standard); Optional: 8mm tube, 1/4-28 UNF bottomsealing threads. Permanent laser-mark identification for 100% traceability. Mass: Dependent on pump size and motor

None of the information supplied by Diener Precision Pumps constitutes a warranty regarding product performance or use. Any information regarding performance or use is only offered as suggestion for investigation for use, based upon Diener Precision Pump's or other customer experience. DPP makes no warranties, expressed or implied, concerning the suitability or fitness of any of its products for any particular purpose. It is the responsibility of the customer to determine that the product is safe, lawful and technically suitable for the intended use. The disclosure of information herein is not a license to operate under, or a recommendation to infringe any patents.

www.dienerprecisionpumps.com


Product Datasheet: Precision Series Piston Pumps

Displacement [milliliters/stroke] vs. Piston Offset Angle					
Offset Angle	Piston Diameter [mm]				
[degrees]	Ø4	Ø6.34	Ø9	Ø12.7	Ø19
0	0.000	0.000	0.000	0.000	0.000
1	0.006	0.015	0.031	0.086	0.208
2	0.012	0.031	0.062	0.172	0.416
3	0.018	0.046	0.092	0.259	0.623
4	0.024	0.061	0.123	0.345	0.831
5	0.030	0.076	0.154	0.431	1.038
6	0.036	0.092	0.184	0.516	1.245
7	0.042	0.107	0.215	0.602	1.451
8	0.049	0.122	0.246	0.688	1.657
9	0.055	0.137	0.276	0.773	1.863
10	0.061	0.152	0.306	0.858	2.068
11	0.067	0.167	0.337	0.943	2.272
12	0.072	0.182	0.367	1.027	2.476
13	0.078	0.197	0.397	1.111	2.679
14	0.084	0.212	0.427	1.195	2.881
15	0.090	0.227	0.457	1.279	3.082
16	0.096	0.241	0.486	1.362	3.282
17	0.102	0.256	0.516	1.444	3.482
18	0.108	0.271	0.545	1.527	3.680
19	0.113	0.285	0.574	1.608	3.877
20	0.119	0.299	0.603	1.690	4.073




The chart at left shows the pump displacement at offset angles varying between 0 and 20 degrees. Select an offset angle and piston size and find the pump stroke displacement (try to select a piston size to deliver the required flow at motor speeds of between 100 and 300 rpm.) Offset angles are defined in the illustration below. The actual displacement may vary depending on pressure, temperature, and fluid type.

The chart below shows the relative displacement of each piston size. The piston size is defined by the diameter of the piston (in mm).

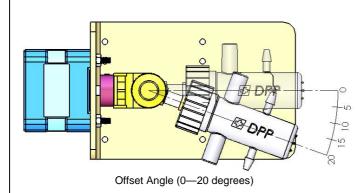

Pump Displacement At Maximum Piston Offset Angle

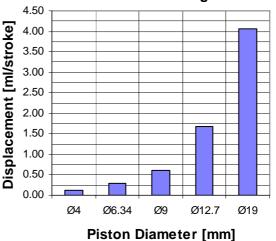

Operating Principle:

